

# SRRF-Stream+ Super-Resolution Microscopy Accessible to All

Fast, reliable & live-cell compatible Super-Resolution

Science has limits imposed by the laws of physics that constrain discoveries and the advance of knowledge. In microscopy, up until the beginning of the 21st century, the diffraction limit of light was an unbreakable barrier. This law of physics imposes that two points could not be resolved (clearly separated) if they are closer together than half the wavelength of light used to view them. In practice, this would mean that light microscopes could only resolve objects/structures that are separated by 200 nm or more. The 200 nm barrier would therefore leave a significant gap of knowledge to be uncovered, since many subcellular structures and organelles are smaller than 200 nm.

In the first decade of the 21st century, the diffraction limit of light was overcome using a number of ingenious microscopy techniques. A new field of discovery was ready to be revealed. Several methods for super-resolution evolved,

allowing imaging beyond the diffraction limit of light. Super-resolution microscopy methods such as STED (stimulated emission depletion microscopy), STORM/PALM (Stochastic optical reconstruction microscopy)/ (Fluorescence photo-activation

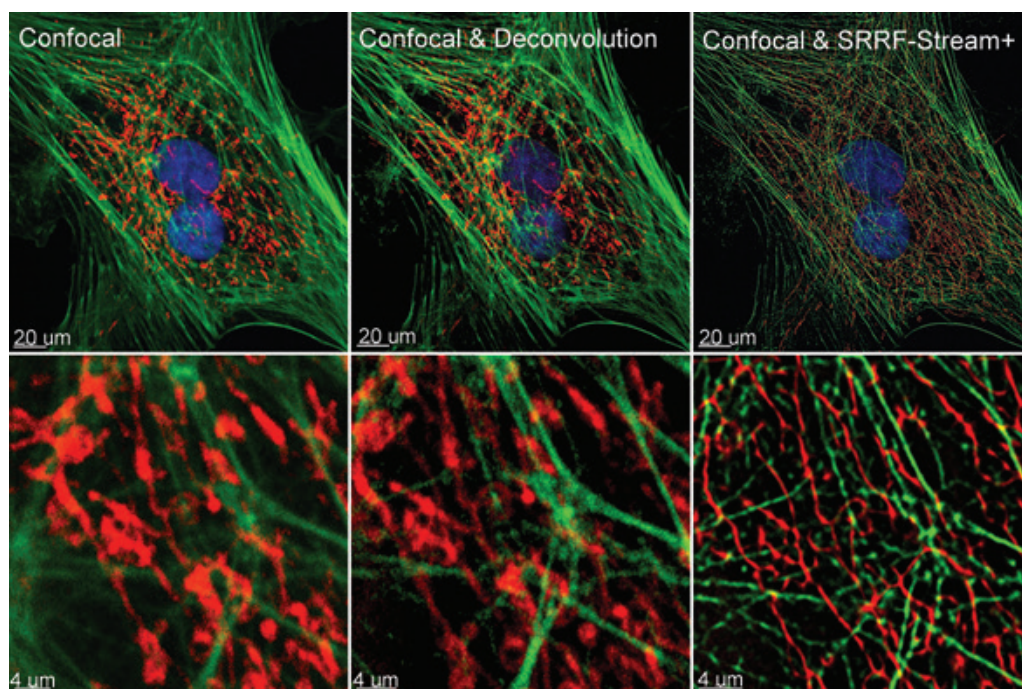


Figure 1 – Comparison between confocal imaging, confocal imaging & deconvolution and confocal imaging & SRRF-stream+. High magnification close ups are present for a better visualisation of the obtained resolution. BPA cells stained with phalloidin, mitotracker and DAPI image with Ixon 888 in confocal mode (no SRRF) and confocal mode with SRRF-stream+.

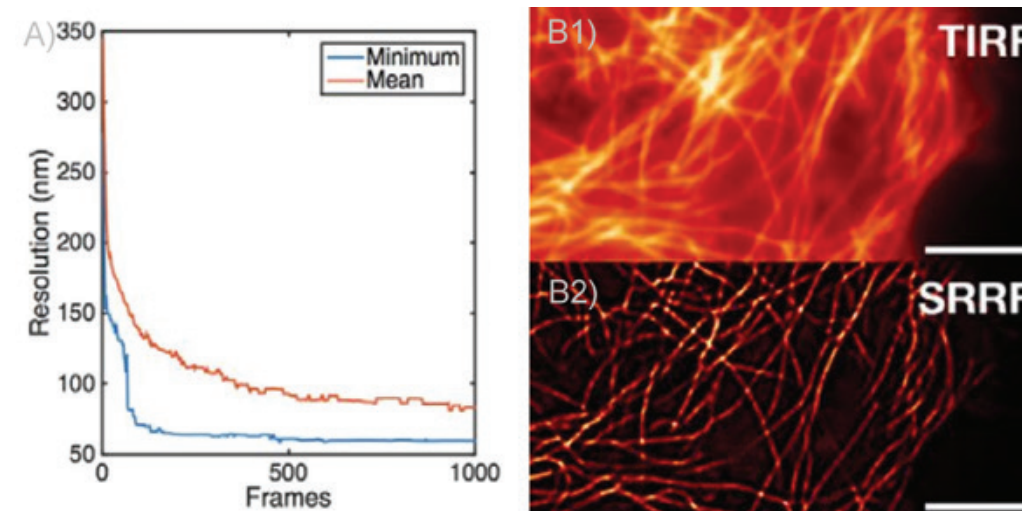


Figure 2- SRRF and SRRF-stream resolution increase with the number of frames acquired (and processed) per time point. A) Graph presenting the resolution increase with the increasing number of acquired frames. As it can be observed there is a high increase up until 100 frames. Mean presents the average of all radial fluctuation correlation analysed (in different Regions of interest - ROI). Minimum shows the maximum resolution achieved for a given ROI. B1 and B2) Actin filaments imaged on Andor Dragonfly using TIRF (B1) and TIRF with on the fly SRRF-stream processing (B2). It can be observed even when acquiring in TIRF modality, that SRRF-Stream provides a significant increase in resolution.

localisation microscopy) and SIM (Structured illumination microscopy) have become available for researchers.

Although the Super-resolution methods have opened a new field of discovery, some limitations on these methods existed. Complex sample preparation, long acquisition times as well as high energy requirements for acquisition, render these techniques inappropriate (if not incompatible) for use with live-cell imaging. Furthermore, the optical requirements and computer power needed render these SR methods extremely expensive and inaccessible to many labs.

More recent advances with spinning disk confocal technology combining optical over-sampling with computational reassignment do achieve a 1.6x improvement beyond the diffraction limit and are suitable for live-cell imaging, but at the significant cost of a reduced field of view, and starting from a lower native resolution. Additionally, it is still inextricably linked to the high-cost hardware.

The Dragonfly spinning disk confocal with its unique combination of pinhole size and custom pinhole

pitch separation has an improved native axial and lateral resolution, delivering better resolution for routine imaging. Then, using photon reassignment (deconvolution) on the acquired images, the final obtained axial and lateral resolution of the dragonfly can be 240 nm and 139 nm respectively.

But what if a researcher needs more resolution? Or does not have a spinning disk confocal microscope? Is there an alternative method that is fast, compatible with live imaging, and that can still acquire super-resolved images at the full field of view and, if required, deep inside the cells? The answer to these questions is yes, with SRRF or SRRF-Stream+ (Figure 1).

In 2016, the Henriques lab developed an alternative approach to super resolution named **SRRF – Super-Resolution Radial Fluctuations** (1). SRRF can be combined with widefield, TIRF or confocal, and the final resolution will depend on the properties of the acquired data sets. Using the SRRF algorithm, the researchers can achieve resolutions in XY up to 50 nm (1). Furthermore, the SRRF algorithm does not require special sample preparation or special fluorophores for acquisition,



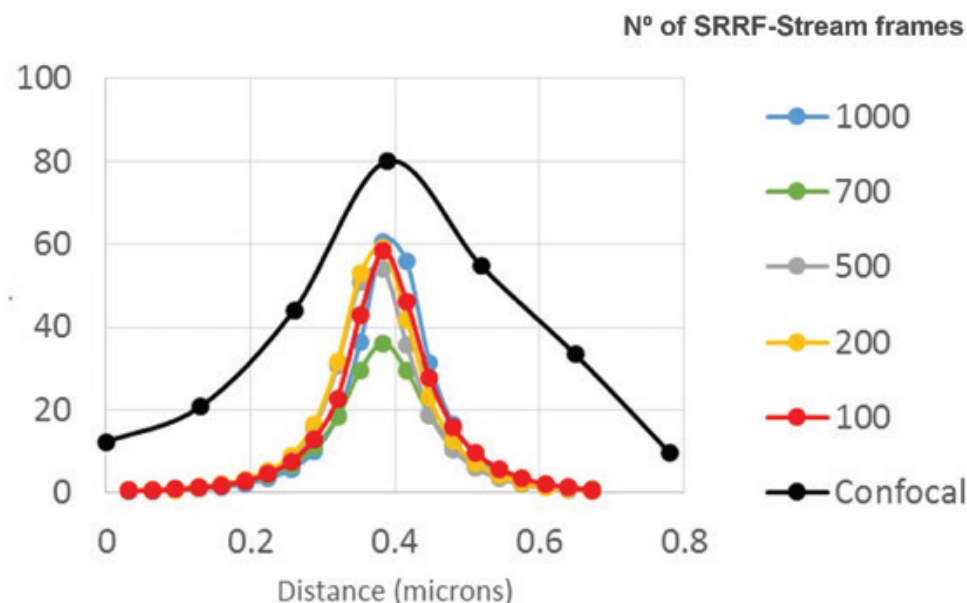


Figure 3- SRRF-stream resolution increase with the number of frames acquired per time point. 100 nm tetraspeck beads were imaged and the FWM was measured. We can observe that with 100 SRRF frames acquired the FWM of the tetraspeck beads is 50 nm. The 100 nm bead can be accurately measured using the SRRF-stream algorithm and acquiring 100 frames. Images were acquired in confocal mode, and in confocal mode with different SRRF-stream acquired image images per point, respectively: 100, 200, 500, 700, 1000. Thanks to Dr Ann Wheeler, Head of the Advanced Imaging Resource at the MRC Institute of Genetics and Molecular Medicine, University of Edinburgh.

being compatible with conventional fluorophores and fluorescent proteins.

Most importantly, a super-resolved SRRF image can be acquired by capturing on average between 20-100 frames (more frames will result in improved resolution) and the energy required for this imaging is in the order of the mW to W per cm<sup>2</sup> range, making SRRF compatible with live-cell imaging. (1, 2) SRRF was initially provided as an ImageJ plugin. In order to obtain a SRRF image, a long acquisition workflow needed to take place and the SRRF-image was obtained by post-acquisition processing of the acquired data in the image J plugin – NanoJ.

In 2018 from collaboration between Andor and professor Henriques, Andor launched **SRRF-Stream**. **SRRF-Stream** is Andor's implementation of the SRRF algorithm that delivers super-resolution images on the fly with the click of a button. As with the original SRRF algorithm, the resolution of **SRRF-Stream** will improve with the number of frames captured per a given time point. This is most striking up to 100 frames,

reaching a more stable increase from 100 frames to 500 frames, and a very modest improvement in resolution from 500 frames onwards (figures 2, 3). Please note that other factors will also affect resolution such as exposure time, Nyquist sampling, radially magnification and the ring radius increase; these should be tested in order to optimise results.

ANDOR's SRRF-Stream was a major improvement allowing live-cell super-resolution, super-resolution deep inside cells or tissues way beyond the top of the coverslip.

SRRF-Stream is available exclusively with **Ixon-EMCCD** cameras, through **micro-manager** and **Fusion** which is the control software for **Dragonfly confocal**. The advantages of SRRF-Stream are (3,4):

- Allows Real-Time super-resolution images, enhancing the workflow and avoids post-processing.
- Breaks the diffraction limit of light by delivering Super-Resolved images with 2 to

6-fold increase in the final resolution (50-150 nm), depending on the imaging mode used (i.e. widefield, confocal or TIRF) and the experimental conditions.

- Use of Low Excitation power Intensities for imaging (mW to W/cm<sup>2</sup>), which results in higher compatibility with extended live cell observations and with minimal impact in cell physiology.
- Compatible with conventional Fluorophore and fluorescent proteins.

### Easy sample preparation.

SRRF-stream is, therefore, a cost-effective solution for any microscope since all of them can be converted into super-resolution microscopes. SRRF-stream is compatible with any imaging modality, namely widefield, TIRF and confocal (Figure 4).

In July 2020 we proudly announced an additional improvement to our SRRF-Stream algorithm: **SRRF-Stream+** (figures 1, 4, 5, 6, 7).

**SRRF-Stream+** goes a step further with increased radiality measurements. On the previous version of SRRF, the radiality measurements were computed in 6 directions and now are computed in 24 directions (Figure 5). This enhancement will improve the result of your super-resolved processing data, and the

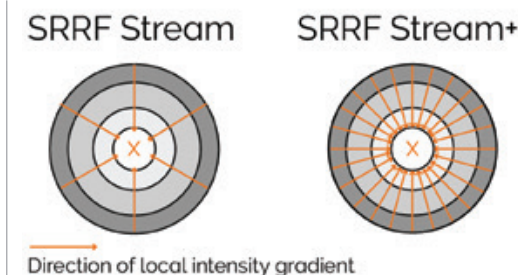


Figure 5 – Radiality computation measurements in SRRF-Stream and SRRF-Stream+. As it can be observed by the image the increased computational radiality measurements will deliver a more accurate result of the SRRF image.

occasional star artefacts previously visible, especially in circular structures such as kinetochores, are now removed (Figure 7).

To get the best from **SRRF-Stream+**, as before, the user will need to acquire images under Nyquist criterium. As previously reported, the major benefit in resolution will be obtained when oversampling by a factor of 2.3 or higher. Still, when oversampling by a factor of 1.5 good results can yet be obtained using the SRRF-Stream+.<sup>1</sup>

When analysing Figures 1, 4, 5 and 7 the benefit on **SRRF-Stream+** is obvious, but one question will come into the mind of many potential users: How is the speed of acquisition affected? If SRRF-stream+ computes radialities in 24 directions instead of 6; is

<sup>1</sup> For more detailed information on how to sample on Ixon cameras please go to Andor original SRRF-stream tech note.

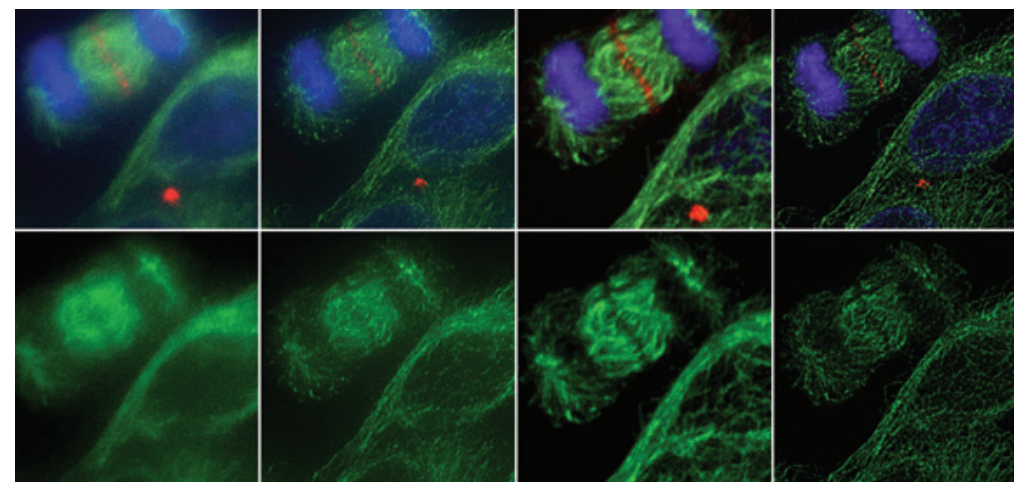


Figure 4 – Comparison between different imaging modes (widefield and Confocal) with and without SRRF-Stream+. Hela cells stained for MLKPI (red), α-tubulin (green-microtubules) and DAPI (blue-DNA) were imaged in the Dragonfly with an Ixon888 camera.



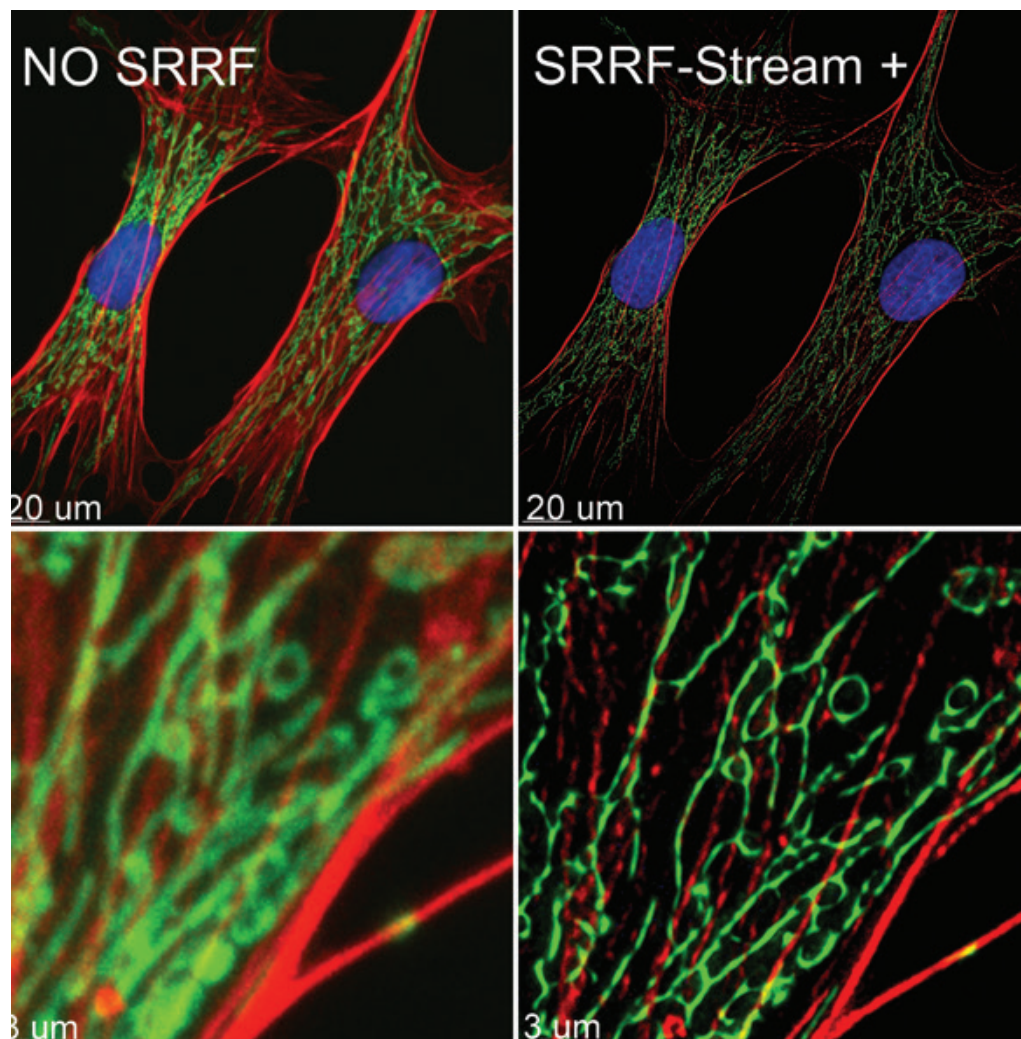


Figure 6 – SRRF-Stream+ delivers High-quality Super-resolved images. BPA cells stained with phalloidin (red), mitotracker (green) and DAPI (blue) where imaged with Ixon 888 Ultra in confocal mode (no SRRF) and confocal mode with SRRF-stream+

SRRF-Stream+ 4X slower than the original version of SRRF? When developing the new SRRF algorithm, Andor took into account that speed of acquisition is also an essential parameter to many researchers. It has been possible to achieve the improved image quality with a minimal impact to the acquisition speeds. To achieve this, CUDA performance has been optimised, perfectly utilising the massive computing power of nVidia graphic card to speed up **SRRF-Stream+** calculations. As a result, a time-lapse of 100 time points with 50 **SRRF-stream+** images per time point (corresponding to 500 images in total) will only vary by 0.5 secs in acquisition when

comparing with our original SRRF-stream<sup>2</sup>.

As for Andor original SRRF-Stream advantages, they are maintained in SRRF-Stream+ but with the improved image quality. Super-resolution imaging processing is still carried out at ultra-fast processing rates (up to 30X faster) than ImageJ SRRF (“NanoJ-SRRF”), and SRRF image acquisition/processing is captured on the fly, in parallel to the acquisition of the data. **SRRF-Stream+** algorithm delivers images with a final resolution between 50-150 nm and due to its low energy requirements is an ideal

<sup>2</sup> Full field of view (1024 X 1024) images acquired with Ixon Ultra

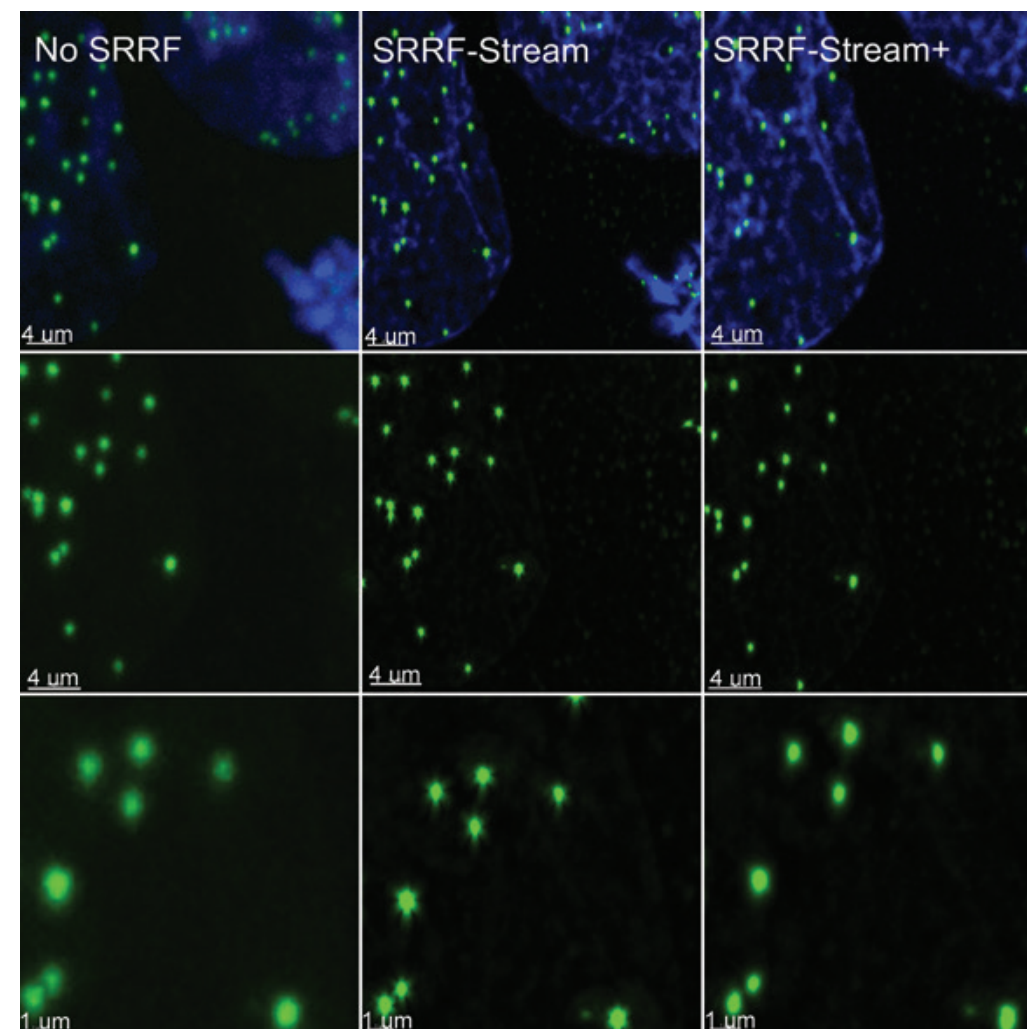


Figure 7 – SRRF stream+ delivers high-quality data. At very high magnifications, in specific structures, such as kinetochores some Star artefact could occasionally appear. SRRF-Stream+ radially computation in 24 directions eliminates the star artefacts. HeLa cells stained for Cenp-A (green-Kinetochore) and DAPI (blue-DNA) were imaged in the Dragonfly with an Ixon888 Ultra camera.

solution for live-cell microscopy. Super-resolved images can be acquired with large fields of view, and importantly super-resolution microscopy is not restricted to cell surface events. With **SRRF-Stream+**, it's possible to acquire a super-resolved image deep inside cell or tissues, and there is no need for special sample preparation.

Since its launch, SRRF-Stream has become widely used for many applications. Examples of using SRRF-Stream in iXon cameras include: to analyse mitochondria trafficking into cell terminus via microtubule destabilisation (5), or analysis of HIF1α

nuclear translocation in mesenchymal stem cells (6). The Dragonfly multimodal confocal system can also be used to acquire SRRF-Stream images, examples are: analysis of cytokinesis using optogenetic tools (7) as well as vesicle trafficking studies on analysing the effects of class II PI3Ks in the regulation of clathrin-dependent pinocytosis (8). Other applications of SRRF-Stream+ include analysis of protein structure at a sub-organelle level, tracking of single molecules inside cells, membrane fusion studies of individual SNARE protein machinery, intracellular skeleton reassembly (changes to actin fibre meshwork).



In conclusion, we have improved our SRRF-Stream algorithm delivering live-cell compatible super-resolution with even better results than before. **SRRF-Stream<sup>+</sup>** delivers super-resolved images of intracellular structures without artefacts, can compensate for the effects of cameras' fixed pattern noise, and delivers high-quality super-resolved images.

Many other applications and discoveries are waiting to be revealed using SRRF-Stream<sup>+</sup>. Do you want to know more about SRRF-Stream<sup>+</sup>? Contact us and we will be delighted to show it in action.

## Bibliography

Gustafsson, N., Culley, S., Ashdown, G. et al. **Fast live-cell conventional fluorophore nanoscopy with ImageJ through super-resolution radial fluctuations.** *Nat Commun* 7, 12471 (2016). <https://doi.org/10.1038/ncomms12471>

Culley S, Tosheva KL, Matos Pereira P, Henriques R. **SRRF: Universal live-cell super-resolution microscopy.** *Int J Biochem Cell Biol*. 101:74-79 (2018). <https://doi.org/10.1016/j.biocel.2018.05.014>

Browne M. **Twelve Reasons Why Your Next Confocal Should Be Dragonfly.** (2017) <https://www.oxinst.com/assets/uploads/downloads/Dragonfly-Twelve-Reasons-Why.pdf>

Coates C, **'SRRF-Stream': Real-Time Super-Resolution in a Camera.** (2017) <https://andor.oxinst.com/assets/uploads/documents/srrf-stream-technical-note.pdf>

Choi, G.E., Oh, J.Y., Lee, H.J. et al. Glucocorticoid-mediated ER-mitochondria contacts reduce **AMPA receptor and mitochondria trafficking into cell terminus via microtubule destabilisation.** *Cell Death Dis* 9, 1137 (2018). <https://doi.org/10.1038/s41419-018-1172-y>

Lee, H.J., Jung, Y.H., Oh, J.Y. et al. **BICD1 mediates HIF1 $\alpha$  nuclear translocation in mesenchymal stem cells during hypoxia adaptation.** *Cell Death Differ* 26, 1716–1734 (2019). <https://doi.org/10.1038/s41418-018-0241-1>

Castillo-Badillo J.A., Bandi A. C., Harlalka S., and Gautam N. **SRRF-Stream Imaging of Optogenetically Controlled Furrow Formation Shows Localised and Coordinated Endocytosis and Exocytosis Mediating Membrane Remodeling.** *ACS Synth. Biol.* 9, 4, 902–919 (2020). <https://doi.org/10.1021/acssynbio.9b00521>

Aung, K.T., Yoshioka, K., Aki, S. et al. **The class II phosphoinositide 3-kinases PI3K-C2 $\alpha$  and PI3K-C2 $\beta$  differentially regulate clathrin-dependent pinocytosis in human vascular endothelial cells.** *J Physiol Sci* 69, 263–280 (2019). <https://doi.org/10.1007/s12576-018-0644-2>

**Claudia Florindo, July 2020**

### Dr Claudia Florindo

*Microscopy Product Specialist, Andor Technology*

Claudia graduated in Applied Chemistry, has a MSc in Human Molecular Biology, a PhD in Cell Biology, and two postdocs in Cell Biology.



With more than 15 years' experience in research, Claudia has worked in several areas like cell biology, HIV studies, oncobiology, endocrinology and neurosciences. In 2010 Claudia joined Universidade do Algarve as a facility manager and three years after Claudia become the director of Advanced Light Microscopy Facility.

In May 2019, Claudia joined Andor Technologies as a Product Specialist in the Microscopy Systems Division.

# Dragonfly

The Most Flexible Imaging Solution Ever!

The game-changer in confocal microscopy. **Andor Dragonfly** multi-modal imaging system delivers instant confocal, widefield, TIRF, dSTORM and SRRF-Stream<sup>+</sup>. all in one system.

The new and improved algorithm SRRF-Stream<sup>+</sup>, computes the radially in 24 directions, further improving the quality of the super-resolved images. Users will benefit from real-time camera-based super-resolution for live-cell imaging across large fields of view, as fast as 1 frame per second (512x512).

SRRF-stream<sup>+</sup> is compatible with conventional fluorophores, with any imaging modality, and even works in thick samples.

